




Simple Perfusion Kit Assembly

8-Channel

ALA Scientific Instruments
60 Marine Street
Farmingdale, NY 11735

T: 631-393-6401 F: 631-393-6407
www.alascience.com

Check before you start:

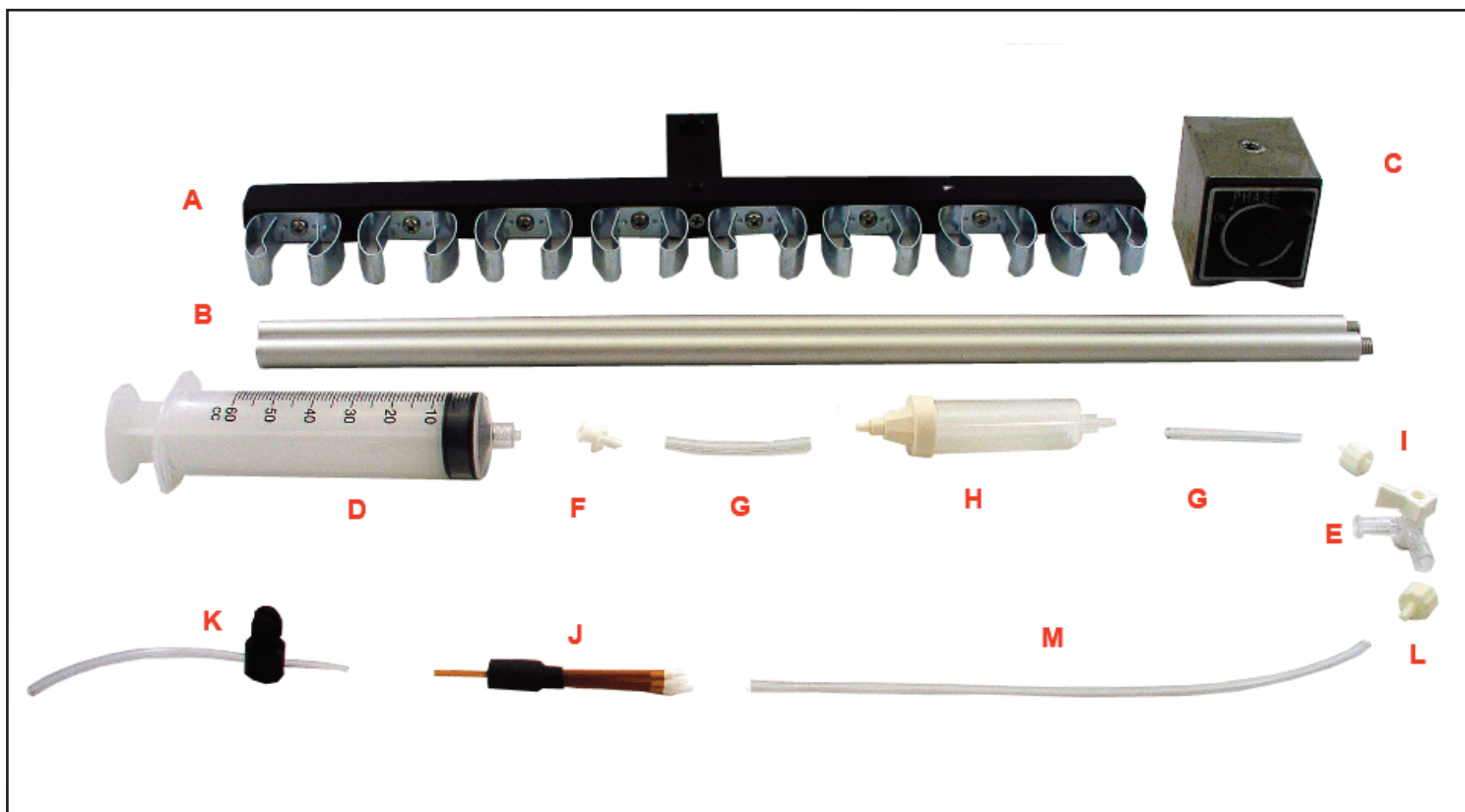

A: Reservoir Bracket x1

B: Magnetic Bar x 2

C: Magnetic Base x 1

D: 60ml Reservoir x 8

E: 3-way Luer Valve x 8

F: Female Barb Luer x 8

G: Connector Tube x 16

H: Drip Chamber x 8

I: Male Barb Luer x 8

J: Minimanifold-8 x 1

K: Control Valve x 1

L: Small Male Barb Luer x 8

M: PVC Tubing x 8

Chamber Assembly:




Connect the chamber, luer valve, luer fittings and tubing as shown above.

Repeat for all 8 channels. Each **M** tube will connect to an input of MMF-8.

Control Valve Attachment and Magnetic Stand:

Attach Flow Control Valve to manifold as shown

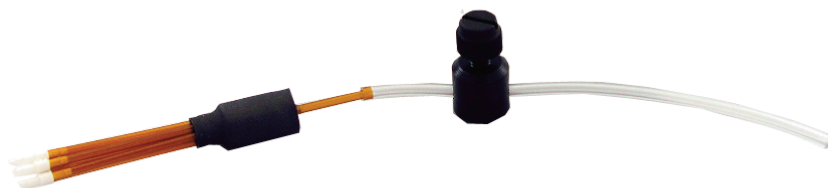

Set up stand as shown in bottom image. Each 60ml reservoir will snap into the clips attached to the bracket.

